




FOR



Up to 1-week Extended Runtime

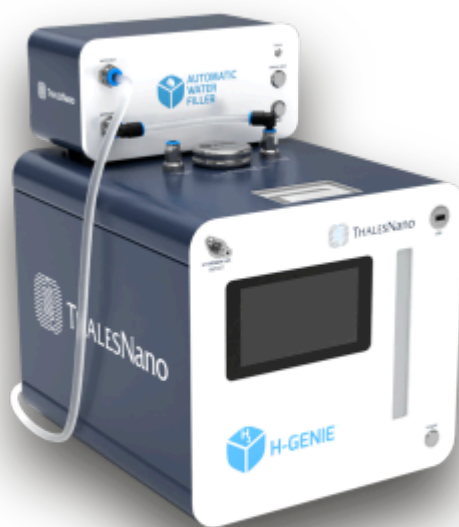

## ENJOY THE SAFETY OF UNLIMITED HYDROGEN GENERATION WITHOUT THE NEED TO MANUALLY MONITOR THE WATER TANK LEVEL

### SAFETY FEATURES

When paired, the **Automatic Water Filler** recognizes when the tank of the H-Genie® starts to run out, and it initiates a **refilling cycle**. Once finished, it **automatically shuts down**. It also monitors the **water quality**, and if the quality drops, it shuts down immediately.

### AUTOMATIC REFILLING ABILITY

- From an external water supply
- Via the built-in water network



“MILLIPORE MILLI-Q” WATER QUALITY REQUIRED



### COMPATIBLE PRODUCTS


- H-Genie® II
- H-Genie® Lite
- THS System Controller®
- THS ReAction® Software



## TECHNICAL PARAMETERS

<b>Recommended environment</b>	Indoor usage
<b>Power requirements</b>	Mains: 100-240V AC, 47-63Hz
<b>Power input</b>	Max. 1500 VA
<b>Unit Dimensions (H x W x D)</b>	150 mm x 305 mm x 187 mm
<b>Weight</b>	5 kg
<b>Outlet parameter</b>	Tubing OD: 10 mm. The output valve can accept any connector with a male thread Press 1/8"
<b>Operating and storage/transport environment</b>	10-35°C, <70% RH

## OPERATION FLOW CHART

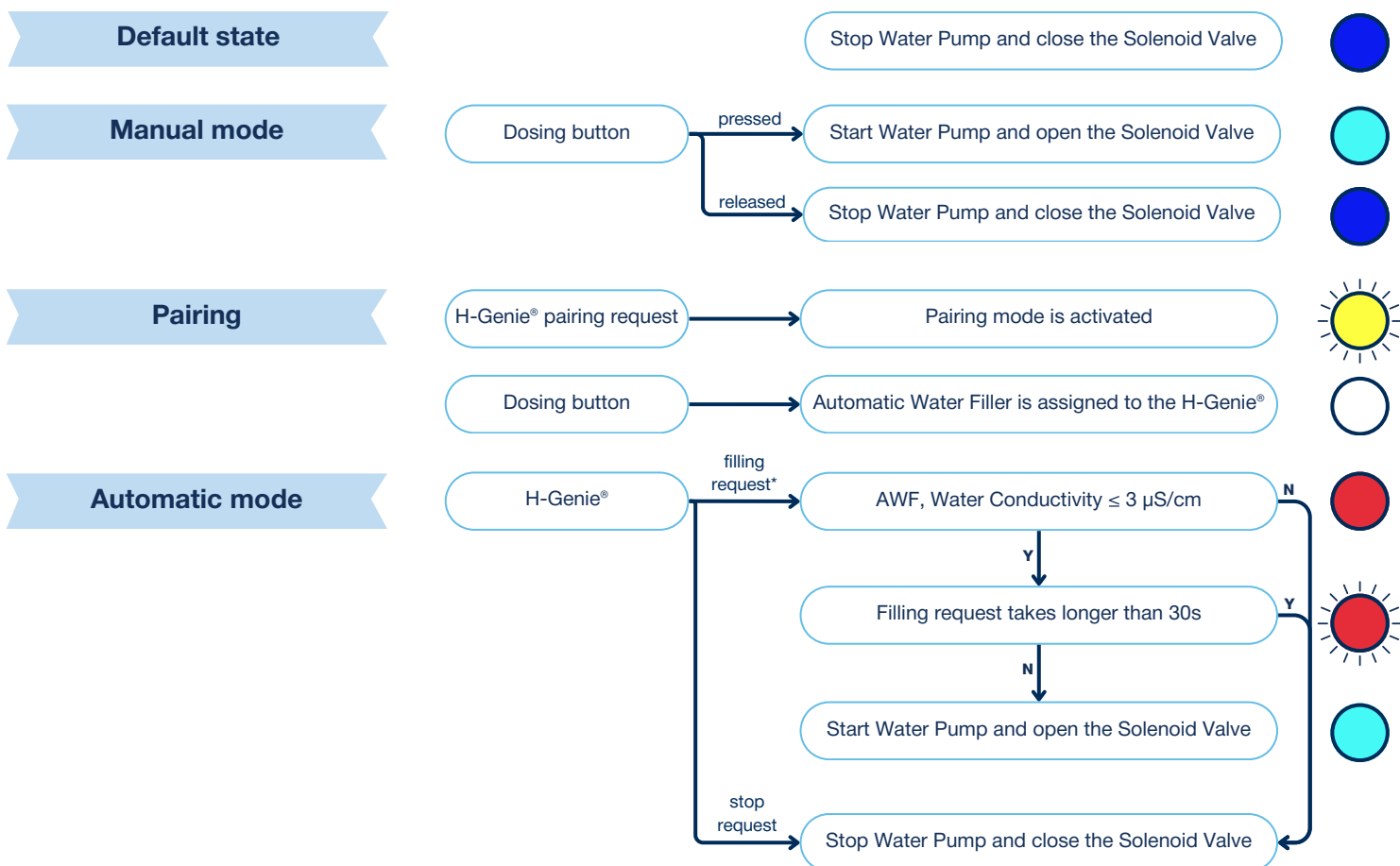

\*the H-Genie® sends water filling request when the upper and the bottom level switches are in down position

For more information, please visit [www.thalesnano.com](http://www.thalesnano.com)



**ThalesNano Inc.**

Zahony str. 7  
 Budapest, ZIP: H-1031, Hungary  
 Email: [sales@thalesnano.com](mailto:sales@thalesnano.com)  
[www.thalesnano.com](http://www.thalesnano.com)

**US Office**

50 S. Penn St. Suite B-2  
 Hatboro PA. 19040, USA  
 Phone: 215-534-3365  
 Email: [USAsales@thalesnano.com](mailto:USAsales@thalesnano.com)

